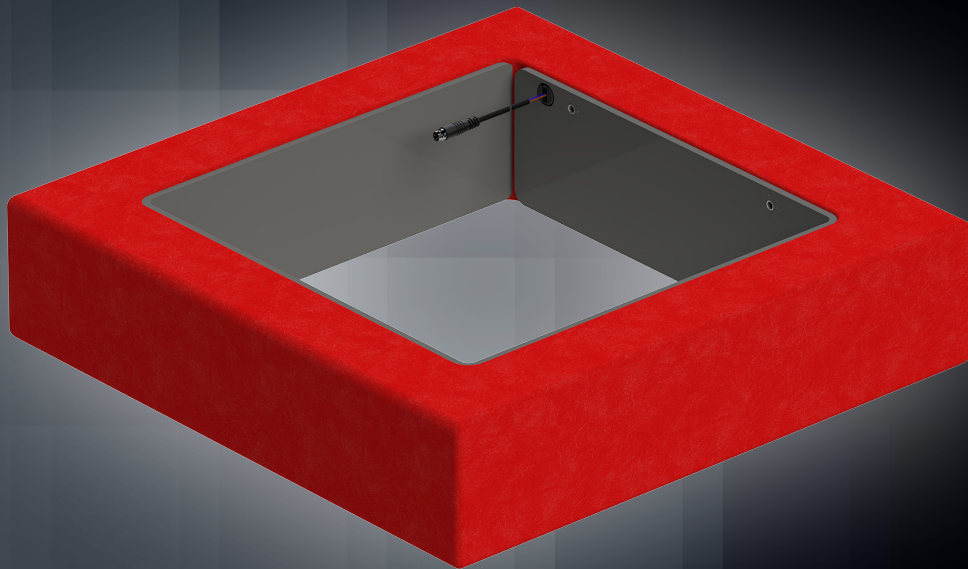
# SENTIR® *bumper*

4 robotics

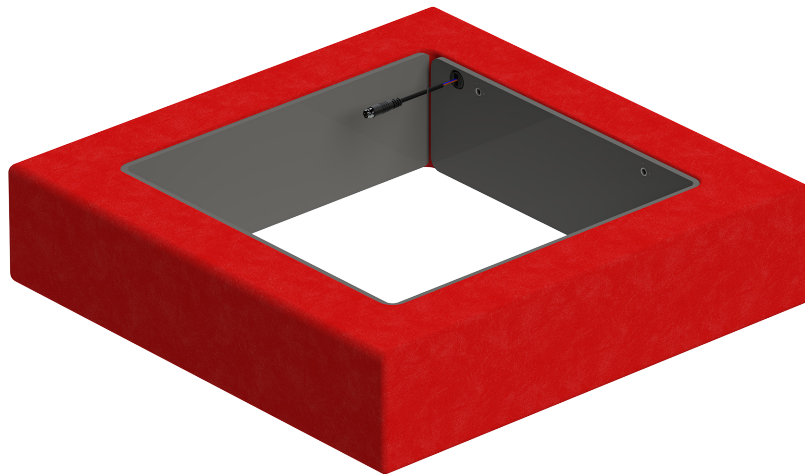
Produkt Information  
Product Information

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## Specifications

SENTIR bumper  
4 robotics

SENTIR bumper 4 robotics

ASO-Sicherheits-  
kontaktpuffer  
ASO-Safety  
contact bumper

für Roboterarme  
for robot arms

**SENTIR**<sup>®</sup>  
*bumper*

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2006/42/EG  
EN ISO 13856-3:2013  
EN ISO 13849-1:2008/  
AC:2009  
2011/65/EU (RoHS)

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## Technical data

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### Mechanical properties

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\*Characteristics for test temperature +20 °C and 100mm/s test speed with Ø80 mm test piece on position C3.

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The response time of the used controller affects the measured actuating distances of the sensor.

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
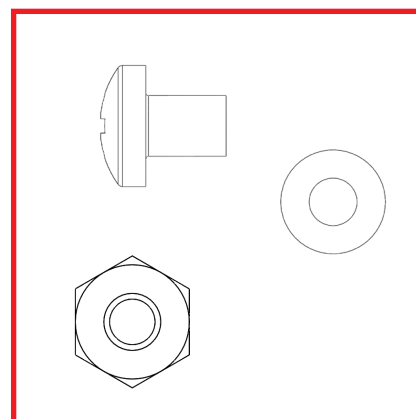
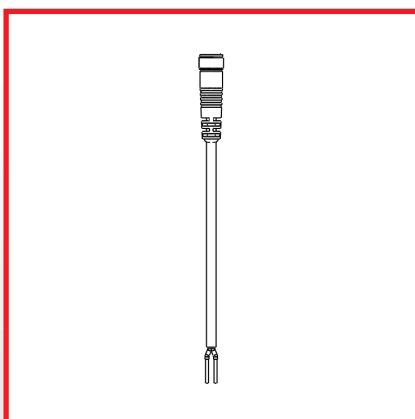
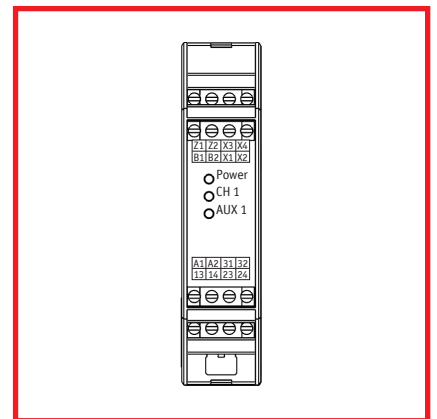
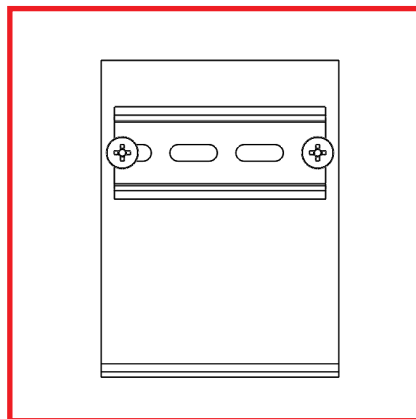
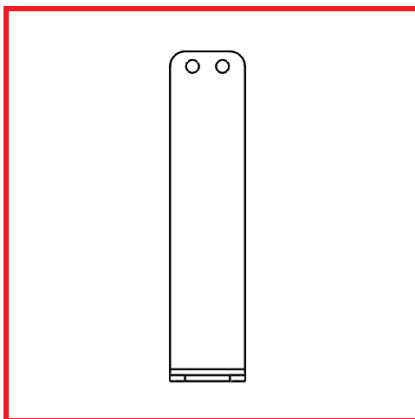
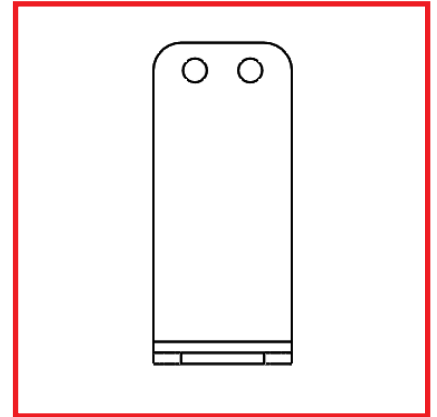
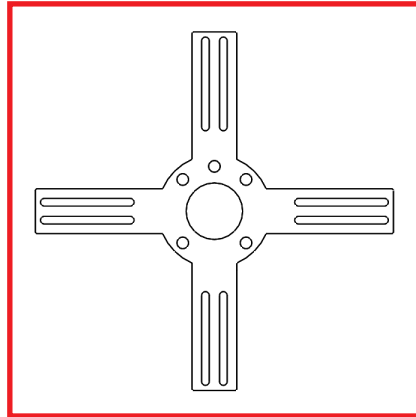
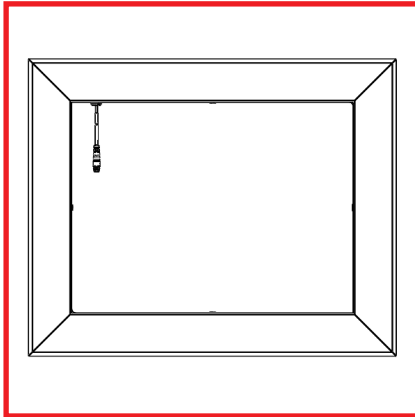
Safety Bumper with mount  
 Extension cable  
 Safety relay with mount  
 Screws

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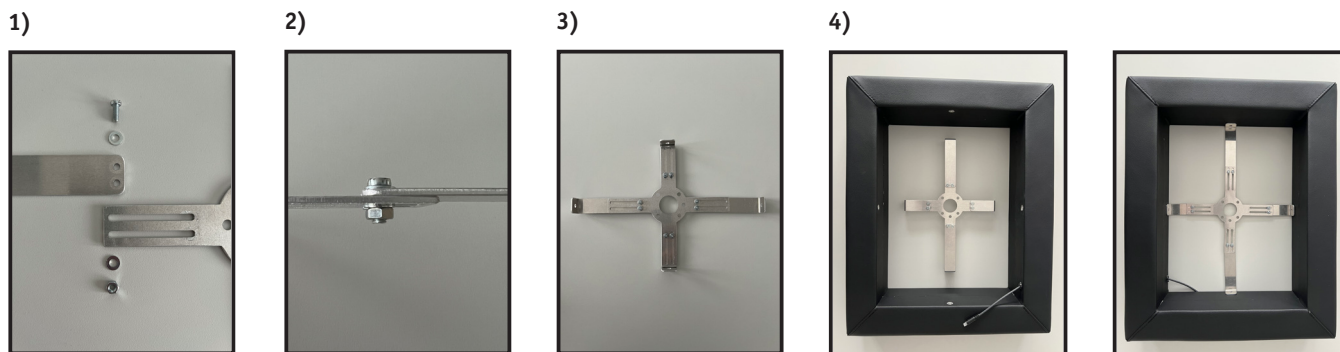Connect the four movable modules to the mounting plate. You will need the standard parts supplied for this. Tighten the standard parts only to 80% so that the holder can be resized.

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Note that the mounting plate is at the top and the movable modules are at the bottom. The screw head lies on the mounting plate with an additional washer (see picture).

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Assembled state of the fixture (picture may differ as there are two different sizes of the movable modules). These are supplied to fit depending on the size of the bumper.

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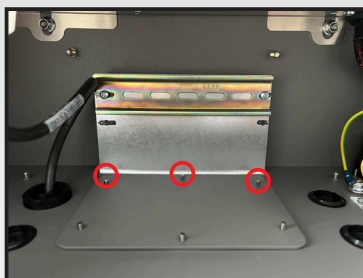
Now pull the movable modules up to the inner surfaces of the bumper. Now tighten the screws 100% tight in the slotted holes. Screw the movable modules into the four threads provided on the bumper using the screws (M4x10) and washers supplied.

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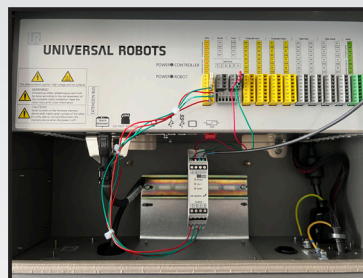
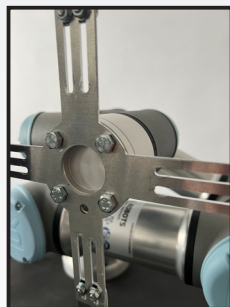
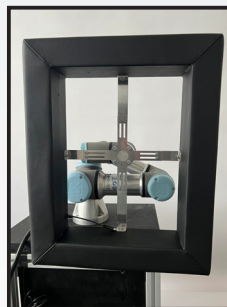
Loosen the six screws and remove the cover plate.

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Slide the ELMON holder under the cover plate onto the screws marked in red as shown in the picture on the left. Make sure that the three holes are on top of each other. Tighten all six screws again.

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Clip the ELMON safety relay onto the DIN rail provided and wire it to the controller as shown on page 8.

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Screw the bumper to the tool holder as shown. Pay attention to the alignment of the centering hole.

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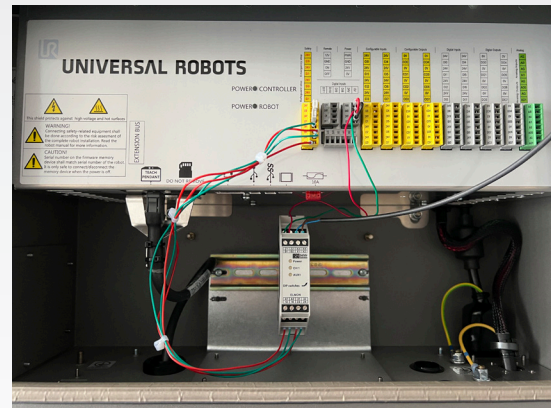
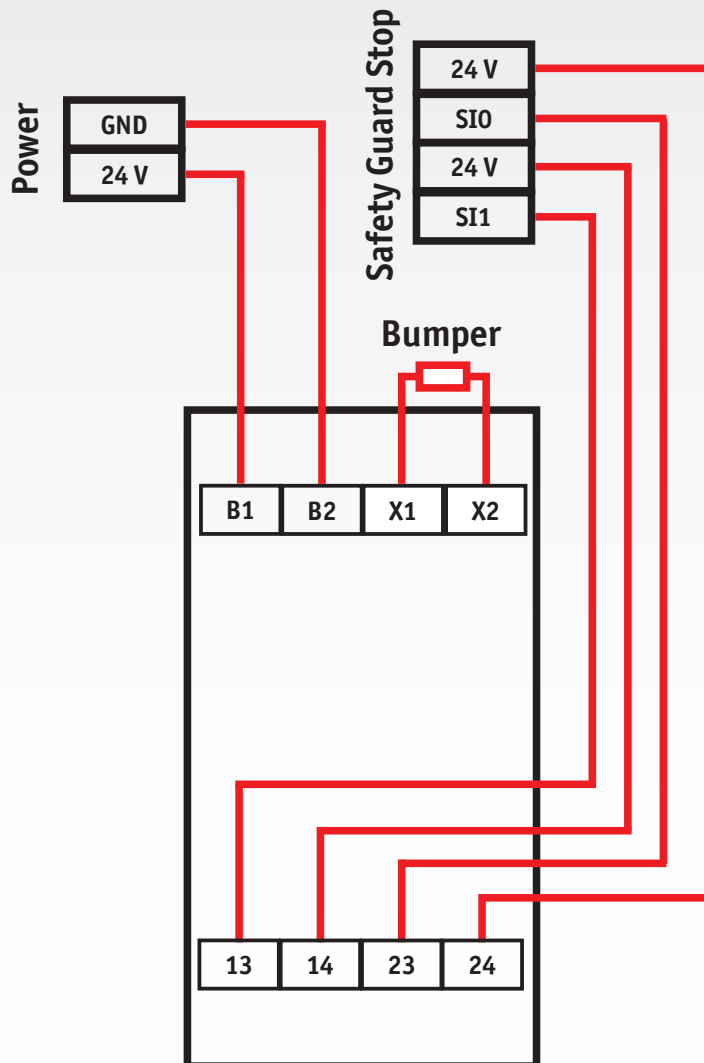
**Before commissioning the robot, please make sure that you adjust the parameters!** Now the bumper can be put into operation.

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Attention, the ELMON relay must of course be set to „Auto Reset“ for the described function in mode 1. Mode 2 must be set to „AUX 1 RL“ off. When the wiring is made as described above, the UR3e control unit can be adjusted as required for the installation. To do this, first go to the „Installation“ group and select the „Safety“ menu item. Via the item I/O, the safety inputs are assigned their functions. In our example we set the function assignment at the line „config\_in[0], config\_in[1]“ to „Reduced mode“. If necessary, we can then parameterize the „Reduced mode“ by going to the „Robot limits“ menu item and adjusting the parameters that are important for us, e.g. the speed.

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We hereby declare that the following products of sensing contact bumpers

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for the combination with safety relays of the model range

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satisfies the relevant essential health and safety requirements of the EC directives and standards listed below on account of its design and construction, as does the version brought to market by us:

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This declaration of conformity does not relieve the designer / manufacturer of the machine from his obligation to ensure that the conformity of the entire machine to which this product is attached satisfies the corresponding EC directive.

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**Manufacturer and attorney of documents**

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#### Headquarters USA

ASO Safety Solutions Inc.  
350 N Frontage Road  
Landing, NJ 07850  
USA

**Phone** +1 973 5869600  
**Fax** +1 973 5861590  
[www.asosafety.com](http://www.asosafety.com)  
[sales-us@asosafety.com](mailto:sales-us@asosafety.com)

#### Headquarters Asia Pacific

ASO Safety Solutions Nanjing Co. Ltd.  
Jinma Lu 3 Maqun Scientific Park  
210049 Nanjing  
CHINA

**Phone** +86 25 856 73990  
[www.asosafety.cn](http://www.asosafety.cn)  
[sales-cn@asosafety.com](mailto:sales-cn@asosafety.com)

DOC0000980 Technical data Rev 00  
as of 31.08.2022  
Technical changes reserved

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ASO GmbH